ESI Electrode Gauge (for Sciex Turbo V Sources)

- cat.# 25832
- $1. \ \ In stall \ electrode \ into \ probe \ per \ in strument \ manufacturer's \ in structions.$
- 2. Lay probe in gauge groove, aligning tip of probe with first line in gauge.
- 3. Use probe adjustment cap to position electrode tip within end of gauge.

Restek, patents and trademarks are the property of Restek Corporation. (See www.restek.com/Patents-Trademarks for full list.)
Other trademarks in Restek literature or on its website are the property of their respective owners. Restek registered trademarks are registered in the U.S. and may also be registered in other countries.

© 2019 Restek Corporation. All rights reserved. Printed in the U.S.A. www.restek.com #735-01-001 Rev. date: 09/19





