ALUGRAM® Xtra

State of the Art **Aluminium Sheets**



200 µm

outstanding wettability easy and reliable excellent separation efficiency



ALUGRAM® Xtra SIL G · aluminium sheets

unmodified standard silica layers on aluminium for TLC

- outstanding wettability for precise colorization results, even with 100 % agueous eluents
- excellent separation efficiency and reproducibilty from lot to lot
- easy and reliable cutting due to an optimized binder system, no flaking of silica

Silica 60, specific surface (BET) \sim 500 m²/g, mean pore size 60 Å, specific pore volume 0.75 mL/g, **particle size 5–17 \mum** Indicator: manganese activated zinc silicate with green fluorescence for short-wave UV (254 nm); special inorganic fluorescence


cent pigment with blue fluorescence for long-wave UV (366 nm)

Binder: highly polymeric product, which is stable in almost all organic solvents and resistant towards aggressive visualization reagents; binder system for ALUGRAM® X*tra* is also completely stable in purely aqueous eluents.

Ordering information

Plate size [cm] Pack of [plates]	2.5 x 7.5 200	4 x 8 50	5 x 7.5 20	5 x 10 50	5 x 20 50	10 x 20 20	20 x 20 25	Thickness of layer	Fluorescent indicator
ALUGRAM® Xtra SIL G · aluminium sheets with standard silica									
SIL G SIL G/UV ₂₅₄	818329	818331	818230.20 818330.20	818261 818360	818232 818332	818362	818233 818333	0.20 mm 0.20 mm	_ UV ₂₅₄

Separation of nutmeg ingredients

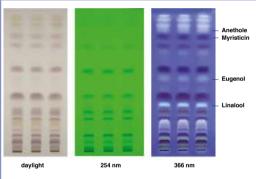

The chromatograms show the following zones with increasing R_t values: linalool (bluish grey), eugenol (yellowish brown), myristicin (reddish brown), and anethole (pink-violet). Other colored zones may appear.

Sample solution: shake 1.0 g freshly powdered drug for 3 min with 4 mL

methanol and filter; apply 10 μL

Layer: ALUGRAM® Xtra SIL G UV₂₅₄
Eluent: toluene – ethyl acetate (95:5, v/v)

Migration distance: 15 cm

Detection: 254 nm: underivatized


daylight and 366 nm: spray with 5% ethanolic sulphuric acid,

1 % vanillic acid and heat to 105 $^{\circ}\text{C}$

MN Appl. No. 403590



Separation of saffron ingredients

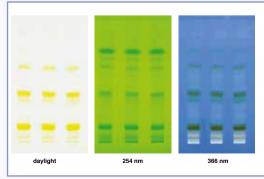

Sample solution: stir 10 mg drug with 50 µL water in a small glass reaction

tube. After 3 minutes add 1 mL methanol and store the solution for 20 minutes in the dark. Afterwards filter through a CHROMAFIL $^{\circ}$ Xtra GF-100/25 filter; apply 10 μ L

Layer: ALUGRAM® Xtra SIL G UV₂₅₄

Eluent: ethyl acetate – 2-propanol – water (65:25:10, v/v/v)

Migration distance: 10 cm

Detection: the sheet is dryed with a hair dryer and analyzed under

daylight, UV 254 nm and 366 nm.

The chromatograms show as main compound Naphtol yellow S.

Other colored zones may appear.

MN Appl. No. 403600

NEW

ALUGRAM® Xtra SILGUR · aluminium sheets standard silica layers with concentrating zone for TLC

- O concentrating zone: valuable aid for manual application and time saving
- excellent separation efficiency
- easy cutting and outstanding wettability


Silica 60, specific surface (BET) ~ 500 m²/g, mean pore size 60 Å, specific pore volume 0.75 mL/g, particle size 5–17 µm

Kieselguhr zone for rapid sample application: because kieselguhr is completely inert towards a large number of compounds, the samples always form a narrow band at the interface of the two adsorbents, irrespective of shape, size or position of the spots in the concentrating zone (see figure). Separation then takes place in the silica layer.

Ordering information

Plate size [cm]	10 x 20	20 x 20	Thickness of layer	Fluorescent indicator		
Pack of [plates]	20	25				
NEW ALUGRAM® Xtra · aluminium sheets standard silica layers with concentrating zone						
SILGUR SILGUR UV ₂₅₄	818412 818422	818413 818423	0.20 mm 0.20 mm	_ UV ₂₅₄		

Save your precious time!

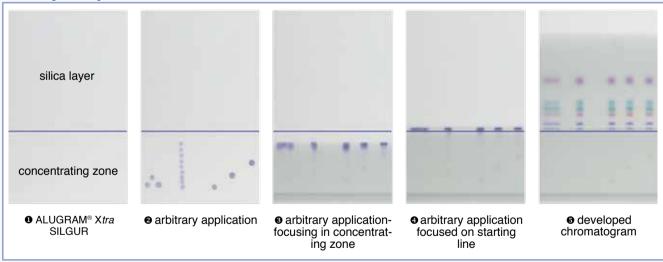

A valuable aid for manual application especially of large volumes of very dilute samples is the concentrating zone •, which consists of a chromatography inactive adsorbent (kieselguhr). The substances to be separated are concentrated to a narrow band in the concentrating zone •. The separation starts at the beginning of the chromatographically active adsorbent silica •.

Concentrating zone as "rapid application zone" - quantitative evaluation of chromatograms is possible, even if samples applied irregular ②. TLC layers with concentrating zone facilitate the handling and also save time in analysis.

Distancia - [---]

ALUGRAM® Xtra Nano-SIL G · aluminium sheets nano silica lavers for HPTLC

- sharper separations in shorter development time and shorter migration distances
- smaller samples and an increased detection sensitivity
- easy cutting and outstanding wettability

Nano silica 60, specific surface (BET) ~ 500 m²/g, mean pore size 60 Å, specific pore volume 0.75 mL/g, particle size 2-10 um

Ordering information

Plate size [cm]	5 x 20	20 x 20	I nickness of layer	Fluorescent indicator
Pack of [plates]	50	25		
NEW ALUGRAM® Xtra Nano-SIL G nano silica layers for HPTLG				
Nano-SIL G Nano-SIL G/UV ₂₅₄	818240 818342	818241 818343	0.20 mm 0.20 mm	_ UV ₂₅₄

NEW

ALUGRAM® Xtra Nano-SILGUR · aluminium sheets

nano silica layers with concentrating zone for HPTLC

- sharper separations in shorter development time and shorter migration distances
- concentrating zone: valuable aid for manual application and time saving
- easy cutting and outstanding wettability

Nano silica 60, specific surface (BET) ~ 500 m²/g, mean pore size 60 Å, specific pore volume 0.75 mL/g. particle size 2-10 µm

Kieselguhr zone for rapid sample application: because kieselguhr is completely inert towards a large number of compounds, the samples always form a narrow band at the interface of the two adsorbents, irrespective of shape, size or position of the spots in the concentrating zone. Separation then takes place in the silica layer.

Ordering information

Plate size [cm]	10 x 10	Thickness of layer	Fluorescent indicator
Pack of [plates]	25		

NEW

ALUGRAM® Xtra · Nano-SILGUR

aluminium sheets nano silica layers with concentrating zone for HPTLC

Nano-SILGUR	818432	0.20 mm	_
Nano-SILGUR UV ₂₅₄	818442	0.20 mm	UV_{254}

www.mn-net.com

MACHEREY-NAGEL



Tel.: +49 24 21 969-0 Fax: +49 24 21 969-199 Fax: E-mail: info@mn-net.com

Switzerland: **MACHEREY-NAGEL AG** Tel.: +41 62 388 55 00 Fax: +41 62 388 55 05

MACHEREY-NAGEL GmbH & Co. KG · Neumann-Neander-Str. 6-8 · 52355 Düren · Germany **MACHEREY-NAGEL EURL** Tel.: +33 388 68 22 68 +33 388 51 76 88 Fax:

MACHEREY-NAGEL Inc. Tel.:

+1 484 821 0984 +1 484 821 1272 Fax: E-mail: sales-us@mn-net.com



KATEN200078/AlugramXtraen3/2/0/06.2013 PD Printed in Germany